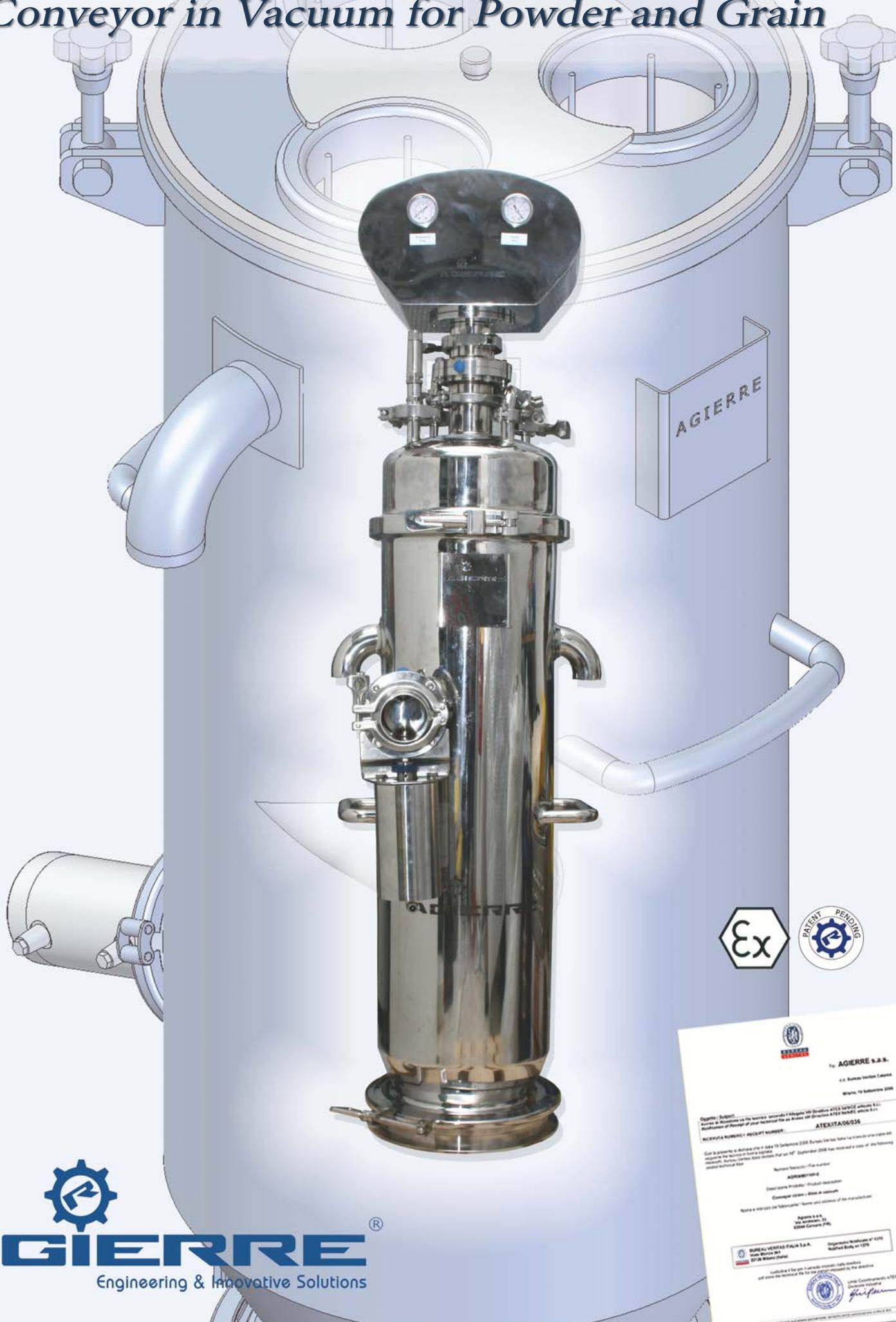
Conveyor in Vacuum per Polveri e Granuli

Conveyor in Vacuum for Powder and Grain

AGIERRE
Engineering & Innovative Solutions

Conveyor in Vacuum per Polveri e Granuli

Powder and Grain Conveyor in Vacuum

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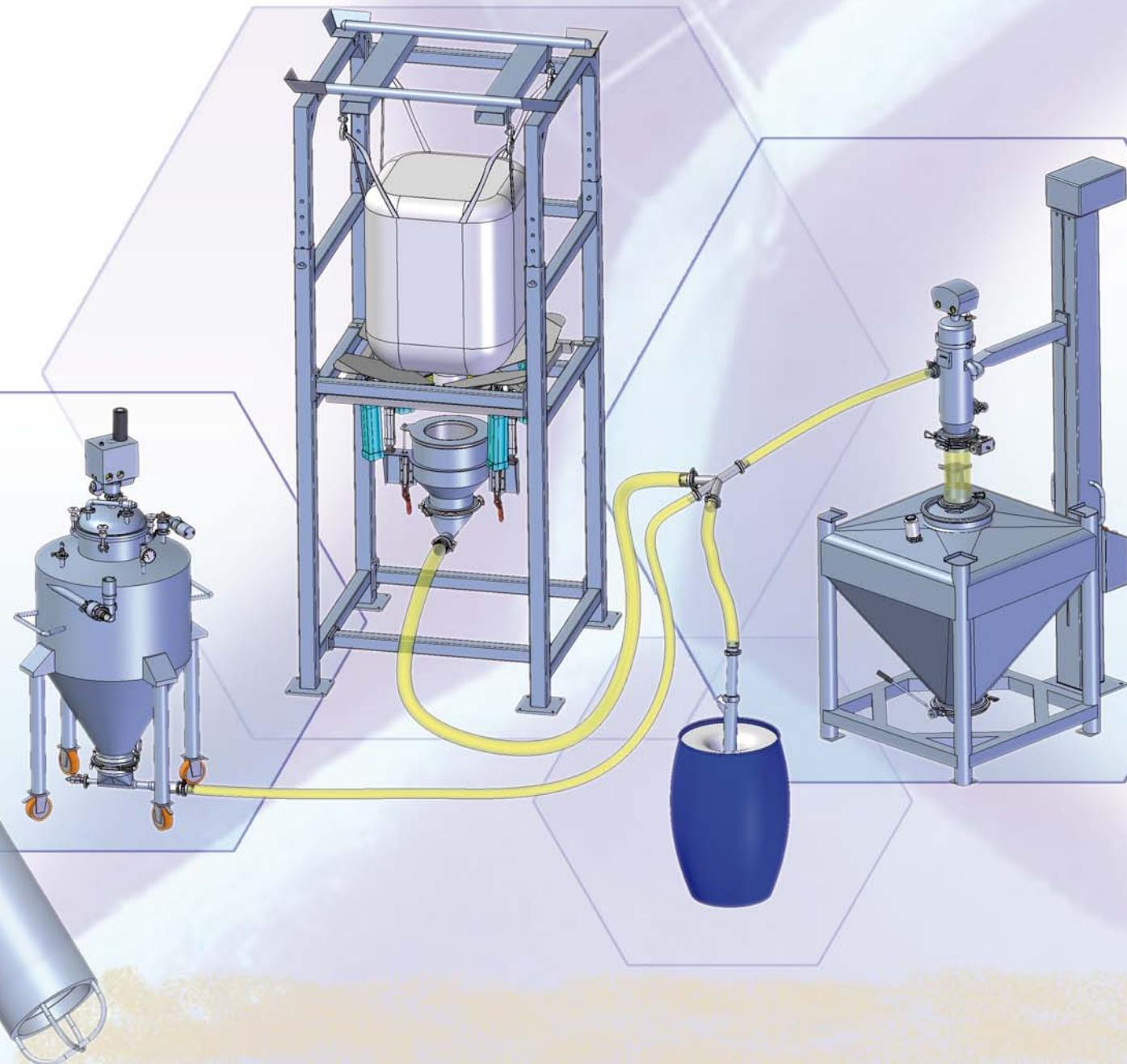
In the last twenty years the experience in this specific field has shown that the Pneumatic Conveyors are the best system in this technology for the handling of powder. Through this system, the powder and grain are transferred from one container to another by means of vacuum aspiration. They are applied both in the productive units (mixers, dryers, fluid bed coolers, blending systems) and in the packing process (packaging machines, table press machines, etc).

A series of injectors, made following the Venturi Principle, generate the vacuum. Compressed air is applied to the injectors generating the depression in the conveyor and consequently, the aspiration of the product. Therefore the product inside it will be separated from the air through a filter. After that, the product falls down thanks to gravity and centrifugal force.

The Vacuum, generated by the pump, allows the handling to a low velocity. Therefore this process guarantees a constantly and consistent handling of the product, without any shock of it. Depending on the product, machines can be built accordingly. The main sectors for the use of the machines are the chemical, pharmaceutical and food industry.

Conveyor in Vacuum per Polveri e Granuli

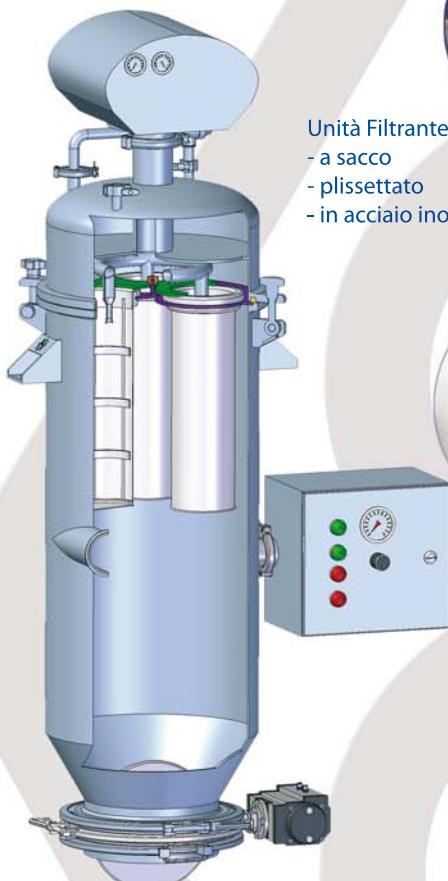
Powder and Grain Conveyor in Vacuum

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This Conveyor has an innovative design for which is deposited a regular patented application. Therefore it is completely modulate concerning the several and different types of connections between pump, cap, machine and the bottom valve. The connections are: flanged (also in the UNI PN16 version). The Conveyor is completely modulate. According to own requirement, these machines can be built accordingly.

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System of Filter:
- bag
- pleat
- stainless steel

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The Conveyor is built following the standard of construction of the medicines of human usage cGMP. This Conveyor has an innovative and peculiar design without any unreachable points or accumulation of product.

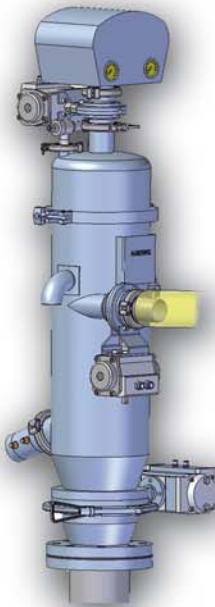
Conveyor in Vacuum per Polveri e Granuli

Powder and Grain Conveyor in Vacuum

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The Conveyor can be used in the applications where there are explosive powders which require during discharge a residual of oxygen inferior of 6%.

In this application there are two pneumatic butterfly valve, one on the incoming of the aspiration and the other on the vacuum line.

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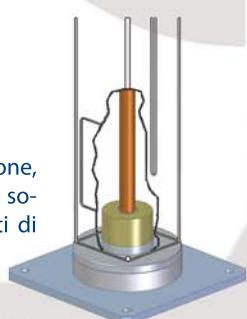
Conveyor with a nitrogen inertization system.

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
This application guar-
tees the stability, avoi-
ding any point of pro-
duct accumulation.

Lifting Column AGR-Lift

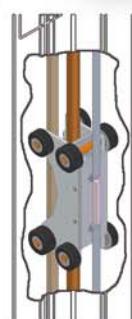The Lifting Column is used in the applications where is necessary to handle the Conveyor. It is in a Stainless Steel AISI 304. This Lifting Column rotates on the main axis of 360° guarantying the right positioning of the lifted object. The mechanical arm has a variable length and a lifting capacity up to 1000 kg. At the inside of the Lifting Column there is the slide which allows, through an alternate movement, the lifting of the mechanical arm. The kinematics coupling "worm/wheel gear" moves the slide. A speed reducer moves the "worm/wheel gear" with a velocity (climb/decline) from 1 to 4 m/min depending on the model.

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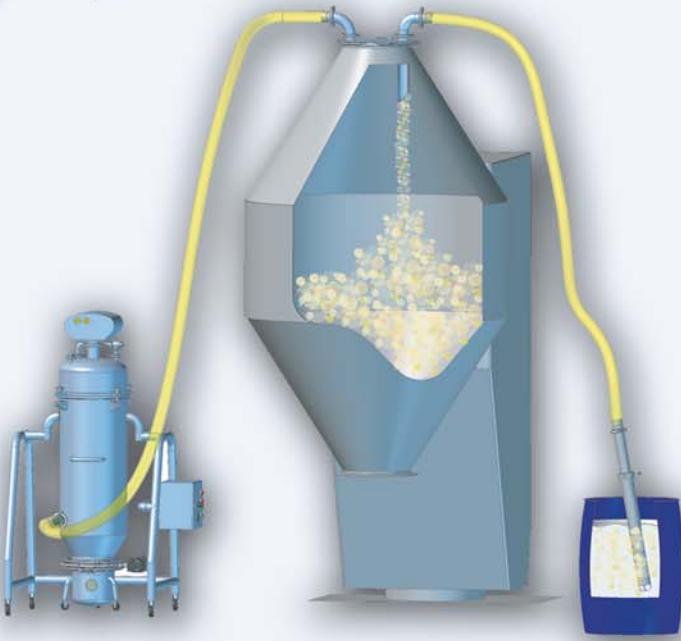
Scarpetta in vacuum *Small aspiration hopper*

Esecuzione Speciale con pompa a vuoto e quadro comandi separati
Application with vacuum pump and control console installed separately from Vacuum Conveyor

**Conveyor in Filter-Vac version for the handling of powder
(in continuous modality)**

The Ciclon-Filter has a sanitary design. Therefore it can be installed directly in this specific area in order to recuperate all residual powder or manually with a container placed directly under the machine or an automatic system with a container in vacuum.

Caratteristiche tecniche / Technical Specifications (Vacuum Conveyor)

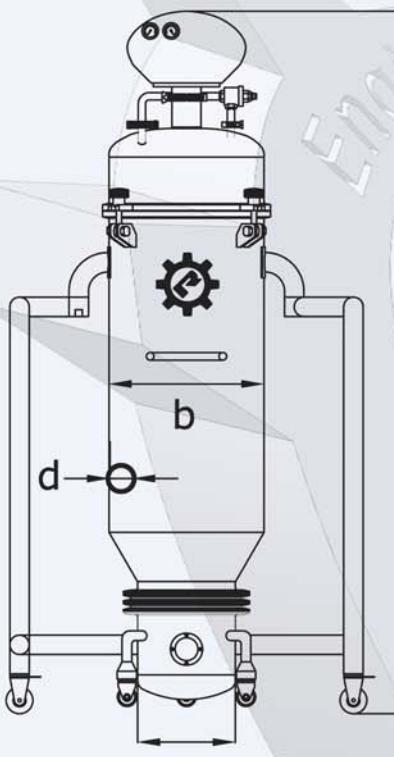
Conveyor in Vacuum per Polveri e Granuli

Powder and Grain Conveyor in Vacuum

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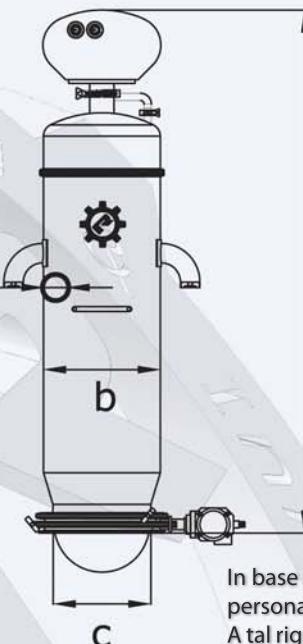
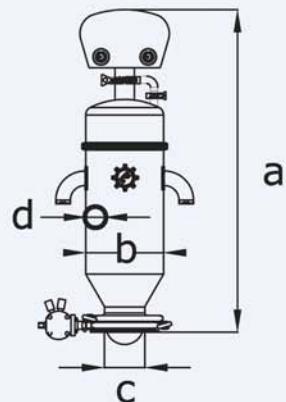
The discharge valve can be realized and customized according to the necessity of the customer.
See the catalogue of our Sanitary Valves.

Dimensioni / Dimension (mm) "Conveyor Agierre"

Tipo / Type		MINIVAC				STARVAC			FILTERVAC		
Modello / Model		AGR150	AGR200M	AGR200	AGR200L	AGR250	AGR300	AGR400	FT250	FT300	FT400
a	Altezza <i>Height</i>	800	840	1070	1070	1320	1410	1570	1180	1680	1840
b	Diametro <i>Diameter</i>	140	200	200	200	250	300	400	250	300	400
c	DN Valvola di Fondo <i>ND Bottom Valve</i>	100	100	150	150	200	250	250	200	250	250
d	Diametro Ingresso <i>Incoming Diameter</i>	1 1/2"	1 1/2"	2"	2 1/2"	2 1/2"	2 1/2"	3"	2 1/2"	2 1/2"	3"

Via Ponte La Pietra, snc
Zona Industriale
I-03043 Cassino (FR) - ITALY
Tel. +39 0776-367 914 - 364 233
Fax +39 0776-364 005
www.agierre.eu - info@agierre.eu